GIMA

MICROSCOPIO
MICROSCOPE
MICROSCOPE
MIKROSKOP
MICROSCOPIO
MICROSCÓPIO
ΜΙΚΡΟΣΚΟΠΙΟ
مجهر

MANUALE D'USO E MANUTENZIONE
USE AND MAINTENANCE BOOK
INSTRUCTIONS DE FONCTIONNEMENT ET ENTRETIEN
BETRIEBS UND WARTUNGS ANWEISUNGEN
MANUAL DE USO Y MANTENIMIENTO
MANUAL DE USO E MANUTENÇÃO
ΕΓΧΕΙΡΙΔΙΟ ΧΡΗΣΗΣ ΚΑΙ ΣΥΝΤΗΡΗΣΗΣ
دليل الإستعمال والرعاية

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ATTENTION: The operators must carefully read and completely understand the present manual before using the product.

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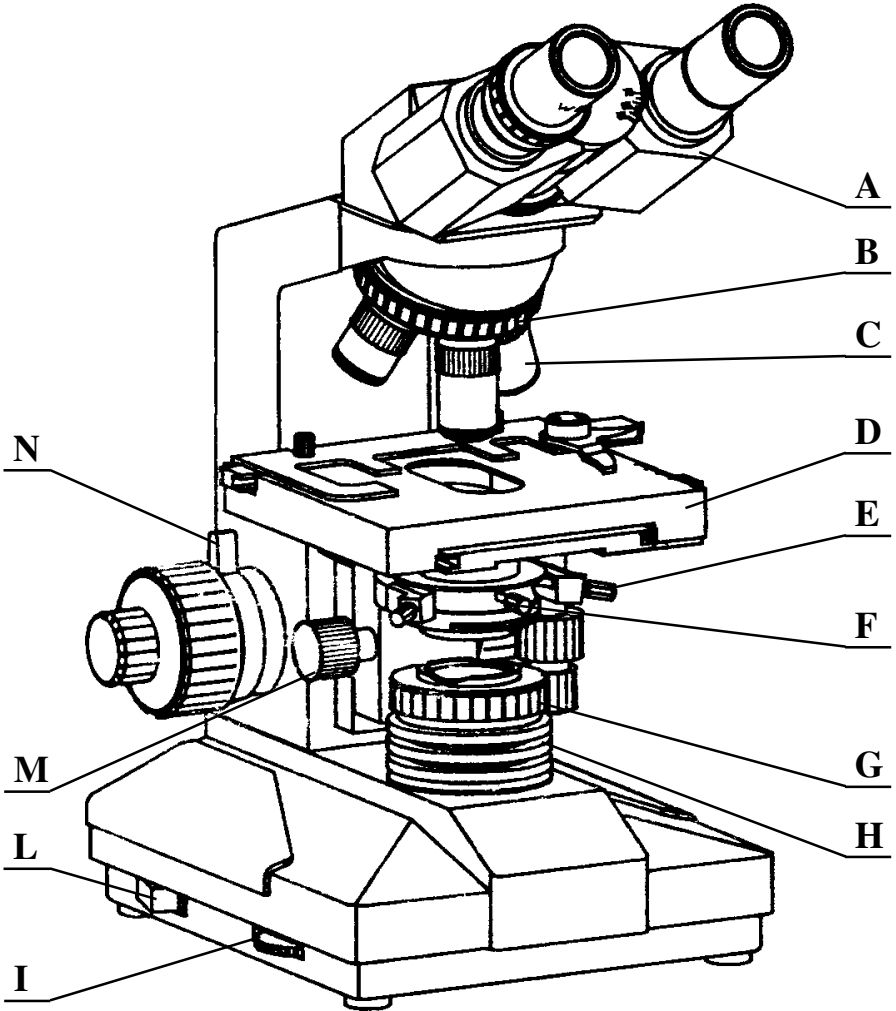
- A Binocular head
- B Nosepiece
- C Objective lenses
- D Plate
- E Condenser adjusting screw
- F Condenser fastening screw
- G Field diaphragm
- H Collector
- I Light adjustment
- L Switch
- M Adjusting knob for condenser height
- N Knob stop

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3. Ingrandimento totale

Obiettivi	4X	10X	40X	100X
Oculari 10X	40X	100X	400X	1000X
Oculari 16X	64X	160X	640X	1600X

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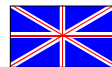
FEATURES

Microscopes GIMA are laboratory equipment supporting the scientific research, for medical use and for didactic use in the schools.

TECHNICAL SPECIFICATIONS

1. Eyepiece lenses

Type	Magnification	Focus (mm)	Field (mm)
Widefield	10X	25	Ø 18
Plan	16X	15.6	Ø 11

2. Objective lenses

Type	Magnification	N.A.	W.D.(mm)	
			Achromatic	Semi-plan
Achromatic or Semi-plan Achromatic	4X	0.1	37.4	23.1
	10X	0.25	6.6	4.1
	40X	0.65	0.64	0.6
	100X (oil)	1.25	0.19	0.38

3. Total magnification

Objective lenses	4X	10X	40X	100X
Eyepiece lenses 10X	40X	100X	400X	1000X
Eyepiece lenses 16X	64X	160X	640X	1600X

4. Condenser numerical opening: NA=1.25;

5. Plate moving field: longitudinal 35 mm, transversal 75 mm;

6. Knob for fine focus: min. increment 0,002 mm;

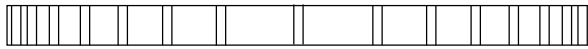
7. Interpupiliary adjustment radius: 53-75 mm;

8. Light source: Halogen lamp with variable intensity from 6V 20W;

9. Power supply: AC 220V 50 Hz or AC 110V 60 Hz;

10. Fungus-resistant

The product, or parts of it, cannot be used for uses different from the one specified in the present manual.

PRESCRIPTIONS

If you move the microscope from a cold to a warm environment, or vice versa, let the instrument acclimate for about half an hour before using it to avoid the formation of condensate.

Do not use the equipment in case it is damaged. Apply to your retailer. Avoid precarious repairs. Repairs shall be carried out with original spare parts only, which shall be installed according to the intended use.

Since the product is made of corrosion-proof materials suitable for the environmental conditions foreseen for its normal use, does not require special care, however it is necessary to store it in a closed place making sure that is protected from dust and dirt to assure its hygienic conditions. Moreover, it is recommended to store the product in a place which can be reached easily by the personnel in case of necessity.

UNPACKING

Always remember that packing elements (paper, cellophane, stitches, adhesive tape, etc.) can cut and/or hurt if they are not carefully handled. They shall be removed with adequate means and shall not be left at the mercy of irresponsible persons; the same is valid for tools used to remove packages (scissors, knives, etc.).

After opening the packages, first of all it is necessary to check all pieces and parts composing the product. Check that they are all present and in perfect conditions.

INSTALLATION

After unpacking the microscope, 180° turn the binocular head:

- Loosen the screw on the binocular head right side
- Gently turn the binocular head by 180° by tightly holding the microscope column
- Tighten the fastening screw again

Each time it is necessary to pack the product, 180° turn the binocular head.

OPERATION

OBSERVATION

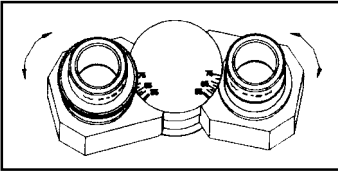
1. Bring the master switch to position "I", to turn the equipment on;
2. Bring the 10X lens to the operating position by means of the nosepiece wheel, then focus the specimen on the plate;
3. Adjust the interpupillary distance and the binocular head dioptré;
4. Adjust the condenser height, the brightness and the diaphragm opening so that to reach a sufficient light effect. When using 4X or 10X lenses, lower the condenser property to reach a symmetrical lighting;
5. To change the objectives, turn the nosepiece wheel and focus again by means of the knob for fine focus. When using the 100X objective lens, place a drop of citron oil between the objective and the specimen.

While turning the nosepiece, pay attention that the specimen does not hit the objectives! The lenses and the specimen can be damaged!


ADJUSTMENT


1. Interpupillary adjustment

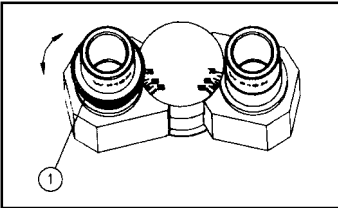Place the specimen on the plate and precisely focus it. Adjust the interpupillary distance until the two views on the right and the left become one.

2. Dioptre adjustment

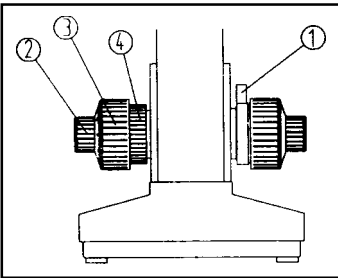
Place the specimen on the plate. Move the 40X lens to the operating position. First look at something with your right eye and focus the right eyepiece by means of the apposite knobs until the image is sharp and clear. Then look at the same thing with your left eye through the left eyepiece and adjust the dioptre 1 until having a clear and sharp image.

3. Rough and fine focus


The instrument makes use of the coaxial focusing device. The tension adjustable knob 4 is used to adjust the tension of the rough focusing knob 3 to prevent the table from lowering due to the gravity. The knob stop 1 is used to avoid that the specimen and the objective can get in contact. The knob 2 adjusts the fine focus.

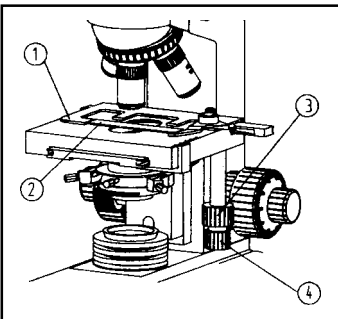While focusing, always pay attention not to touch the specimen with the objective lens!

4. Plate

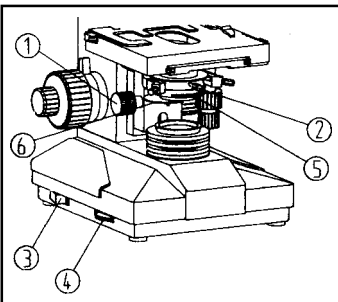The functional holder for the specimen 1 on the plate is used to fix the glass slides 2, the knobs 3 and 4 are coaxial and are used to move the plate along the horizontal axis.


5. Moving condenser

The condenser can be lifted or lowered by turning the apposite knob 1. The condenser can be easily removed by loosening the fastening screw 2. The filters can be fixed to the apposite holder.

6. Master switch and brightness adjustment


Turn the master switch 3 on and adjust the brightness by means of the apposite wheel 4 until the image is clearly visible.

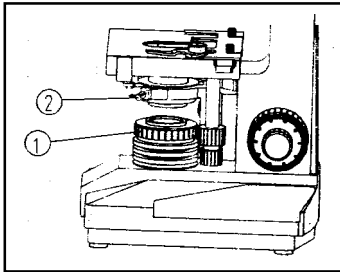Do not leave the brightness wheel to maximum for a long time to avoid reducing the lamp lifetime.

7. Diaphragm field adjustment

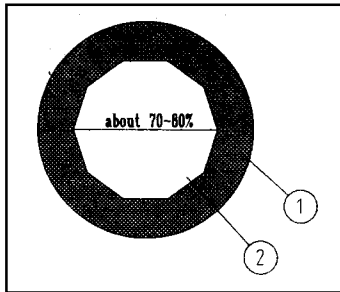Turn the instrument on, place the specimen on the plate, bring the 10X objective lens to the operating position, and watch with 10X eyepieces. Turn the knob to adjust the condenser height to have the diaphragm field image. Focus the diaphragm and the optical axes by means of the adjusting screw **2**. Rotate the diaphragm ring **1** when the diaphragm field is superior to the eyepieces one. Follow the same procedure for the 4X objective lens.

8. Diaphragm opening

The diaphragm opening lever can be rotated to adjust the numerical opening of the lighting system. Remove the eyepieces and watch through the tube. If the diaphragm image is off-centre with reference to the objective **1**, adjust it by means of the adjusting screw. Regulate the diaphragm opening to obtain an image with a good contrast. Usually adjusting the opening diameter of the diaphragm image **2** at 70-80% of the objective it is possible to reach a good result. As a general rule, the wider the diaphragm is opened, the higher the image brightness and resolution; the narrower, the higher the field depth and contrast.

MAINTENANCE

The microscope GIMA is a precision optical instrument and must be handled with great care:

- Place the microscope in dry and clean environment, avoiding sudden changes in temperature.
- When the microscope is not used, always protect it with a dust-proof cover. If you think you do not use it for a long time, replace it into the apposite case supplied with the microscope. For the best preservation of the objective lens and eyepieces, it is advisable to use dehumidification agents.
- In case you need to move the microscope, hold it by the holder column. Do not seize it by the eyepiece or objective lenses.

Before cleaning the microscope, unplug it from power supply!

Never clean the optical surfaces with paper tissues and cloths!

Do not use pure alcohol to clean optical components!

Do not dismantle and never clean the internal surfaces of the microscope optical components!

Remove the oil from the 100X lens after each employment!

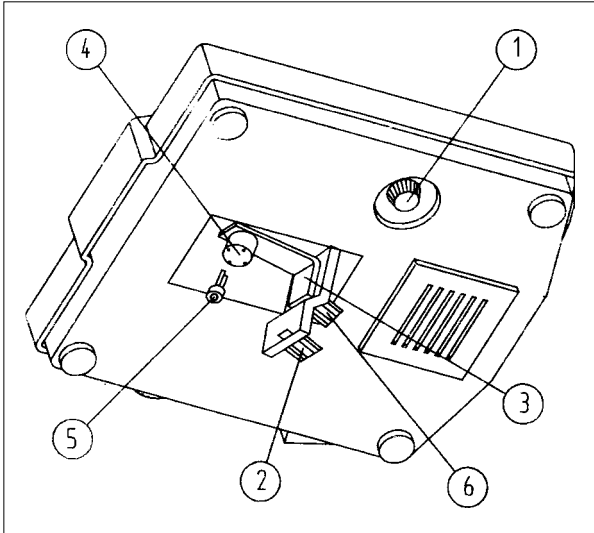
The microscope metallic parts can be cleaned with a slightly wet cloth.

The optical components (objective lenses, eyepieces, condenser) must be daily cleaned as follows: blow away the dust by means of a little pump or compressed air bomb. Alternatively it is possible to remove the dust with a clean brush with non-abrasive bristles.

To thoroughly clean the optical components (objective lenses, eyepieces, condenser) and to remove the oil from the 100X objective, make use of a wad of cotton-wool soaked with a alcohol/ether solution or diethyl benzene.

LAMP REPLACEMENT

1. Turn the microscope off and unplug it.
2. Remove the eyepieces and put them in their case.
3. Incline the microscope, loosen the screw **2** of the cover **3** in the middle of the microscope bottom and remove it.
3. Draw the old lamp out of its seat **4**.
4. Insert the new lamp **5** into its seat **4** and check that it is correctly inserted.
5. Clean the new lamp with alcohol.
6. Replace the cover **3** on the bottom and fasten it by means of the screw **2**.
7. Plug in and turn on the microscope, place the objective and the lenses, adjust the condenser height, and make the light enter the visual field. If the spot is off-centre with reference to the visual field, loosen the screw **6** and

move the lamp base **4**, until it is in the centre, then tighten the screw **6**.

FUSE REPLACEMENT

1. Unscrew the fuse-holder **1**.
2. Draw the old fuse out of the fuse-holder and replace it with a new one.
3. Screw down the fuse-holder **1**.

Fuse specifications: Diameter 5 mm. - 0.5A

Only use fuses compliant with the specifications! Different values could damage the product!

110V POWER ADJUSTMENT

Before carrying out any operation, disconnect the appliance from the mains and wait at least 15 minutes to allow the capacitors to discharge. If you have doubts on the actual network voltage do not perform adjustments! A wrong voltage setting can cause device failure and risk of fire or electric shock!

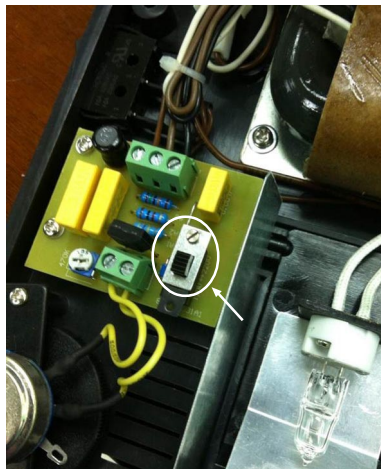
The microscope supplied is set by default to the 220V mains power supply. There is a switch inside the base of the microscope to adjust the voltage supply to 110V.

To adjust, open the microscope base by unscrewing the 4 screws in the corners and move the switch shown in the picture.

If you then need to use the microscope with 220V voltage, remember to move back the switch before connecting the device to the mains!

Disposal: *The product must not be disposed of along with other domestic waste. The users must dispose of this equipment by bringing it to a specific recycling point for electric and electronic equipment.*

For further information on recycling points contact the local authorities, the local recycling center or the shop where the product was purchased. If the equipment is not disposed of correctly, fines or penalties may be applied in accordance with the national legislation and regulations.

GIMA WARRANTY CONDITIONS

Congratulations for purchasing a GIMA product.

This product meets high qualitative standards both as regards the material and the production. The warranty is valid for 12 months from the date of supply of GIMA.

During the period of validity of the warranty, GIMA will repair and/or replace free of charge all the defected parts due to production reasons. Labor costs and personnel traveling expenses and packaging not included.

All components subject to wear are not included in the warranty.

The repair or replacement performed during the warranty period shall not extend the warranty.

The warranty is void in the following cases: repairs performed by unauthorized personnel or with non-original spare parts, defects caused by negligence or incorrect use.

GIMA cannot be held responsible for malfunctioning on electronic devices or software due to outside agents such as: voltage changes, electro-magnetic fields, radio interferences, etc.

The warranty is void if the above regulations are not observed and if the serial code (if available) has been removed, cancelled or changed.

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